
Eyewash bottles BARIKOS

First Aid Measure !




Eyewash bottles are used as a first aid measure to rinse the eyes in order to remove penetrated substances, e.g. dust or chemicals. They should always be at hand and ready for use at hazardous workplaces.

The types BARIKOS and BARIKOS K are designed to be filled with clear water.

The BARIKOS KS and Mini BARIKOS KS models are maintenance-free versions of eyewash bottles. Containing gamma-radiation-treated water, they can be kept ready for use for at least two years in the sealed state. Both models are ideal for keeping in first aid boxes and sets. BARIKOS K and KS eyewash bottles comply with the standard DIN 12930.

8-450 EN 05/2005

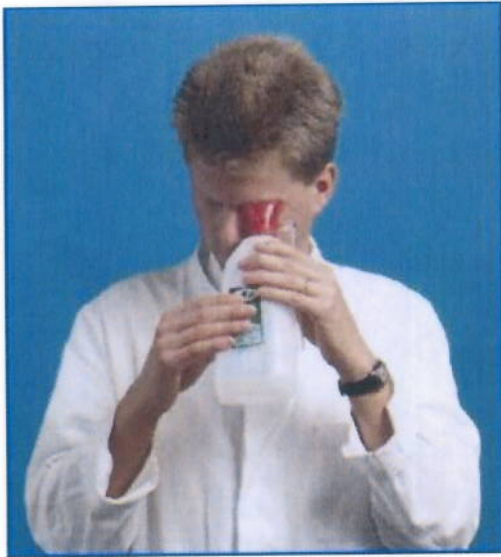

Using an eyewash bottle in the upright position we recommend the type BARIKOS (vol. 620 ml). The user only is obliged to turn his head a little bit downwards coming into good contact to the spray head of the bottle.

BARIKOS: P/N: 954000

The Mini BARIKOS eyewash bottle (vol. 175 ml) is available for persons, being frequently present in different working areas. It can be easily carried along in your workdress pocket.

Mini BARIKOS: P/N: 954003

The BARIKOS K and KS eyewash bottles (vol. 620 ml) can be used immediately even when the injured person is lying down.

As soon as the bottle is tilted downwards, a patented sliding valve opens and the washing fluid reaches the spray head unhindered.

BARIKOS K: P/N: 954002

A protection cover tightens the spray head of the eyewash bottles BARIKOS KS in a hygienically way and without any leaking. The unbroken seal guarantees proper function. Shelf life of the eyewash bottle always is printed on a sticker on front- and backside of the bottle.

BARIKOS KS: P/N: 954004

Mini BARIKOS KS: P/N: 954005



Keeping eyewash bottles everytime in the "ready for use" stage, we recommend the following possibilities:

- fixation support, made of plastic coated steel wire, designed for one eyewash bottle. P/N 154808
- plastic cabinet for wall assembly for one eyewash bottle; the transparent cover is marked with the First-Aid-Symbol. P/N 154812
- plastic cabinet for wall assembly for two eyewash bottles; the transparent cover is marked with the First-Aid-Symbol. P/N 154809

Data Sheet: Eye wash bottle BARIKOS KS

BR-Article-No.:	954004
Trade name:	BARIKOS KS
Product-description:	Eye wash bottle according EN 15154-4:2009-07 with filling free of germs
Kind of product:	First-aid equipment for eye rinsing
Kind of use:	Eye rinsing in standing up-, sitting- and lying position as first aid self help measure to wash out alien elements or substances
Description:	Transparent plastic bottle with printed on pictogram, with red spray head and white cap, internal ascending tube with slide valve, drain tube outside, protection seal
Contents:	Drinking water (ph-value: 7 – neutral), 620 ml, free of active germs by treatment with gamma rays Limited good for use period: 2 years, when seal is unbroken only
Materials:	Bottle, protection cap: Polyethylene Tubes (inside and outside): Polyethylene Spray head, angle piece, slide valve body: Polystyrene Slide valve ring: Polymethylene oxide Washer: Nitrile rubber
Measurements:	approx. 90 x 70 x 250 mm
Weight:	approx. 750 grams
Marking:	BARIKOS KS – EN 15154-4 - 620 ml – Charge „xx“ (Charge-No.) – „yyyy/mm“ (end of “good for use“-time) – 954004 (Article-No.)
Handling/Storage:	At room temperature! Protect from cold, heat and direct solar radiation!
Warning advice:	Eye rinsing does not replace an ophthalmological examination

S 8-501 20141201